

WizScript™ III Reverse Transcriptase

RUO For Research Use Only

REF W2303

DESCRIPTION

WizScript™ III Reverse Transcriptase is a genetically modified M-MLV Reverse Transcriptase which exhibits RNA or DNA dependent DNA polymerase, but lacks ribonuclease H activity. This enzyme can synthesize a complementary DNA strand initiating from a primer using RNA or DNA templates. Removal of the RNase H activity results in an increase of full-length cDNA products. The RNase H activity degrades the RNA part of the cDNA/RNA hybrid following reverse transcription, which may prevent its inhibition during subsequent PCR steps.

KIT CONTENTS

Contents	W2303	W2303-5
WizScript™ III Reverse Transcriptase	10,000 U	50,000 U
5X Reaction Buffer	500 µl	2,500 µl

STORAGE BUFFER

50 mM Tris-HCl, 150 mM NaCl, 0.1 mM EDTA, 1 mM dithiothreitol, 0.1% NP-40 Alternative, 50% glycerol, pH 7.6 (25°C).

5X REACTION BUFFER

250 mM Tris-HCl, 375 mM KCl, 15 mM MgCl₂, 50 mM dithiothreitol, pH 8.3 (25°C).

UNIT DEFINITION

1 unit is defined as the amount of enzyme required to incorporate 1 nmol of dTTP into acid insoluble material in 10 minutes at 37°C using poly r(A)/oligo (dT) as a substrate.

STORAGE CONDITIONS

• Store all components at -20°C in a non-frost-free freezer.

PROTOCOL

1. mix 1-5 µg RNA with
 - 1 µl of 100pmol/µl oligo-dT₁₅₋₂₅ or random primers,
 - or 1 µl of 10pmol/µl gene-specific primer per µg of RNA.
2. Incubate the mixture at 70 °C for 5-10 min
3. Cooling on ice
4. Add the following components in the order listed to a PCR tube

Components	Volume	Final conc.
WizScript™ III Reverse Transcriptase	1 µl	10 U
5X Reaction buffer	4 µl	1 X
2.5 mM dNTP mix	4 µl	200-500 µM
RNase inhibitor (40 U/µl)	0.5 µl	* optional
ddH ₂ O	Up to 20 µl	

5. Mix well and spin briefly.
6. Incubate at 42-50°C for 30-60 min.
7. Incubate at 95°C for 5 min. for reaction stop.
8. Store products at -20°C or proceed to next step.

NUCLEASE CONTAMINATION TESTS:

Single-Stranded Exonuclease Activity

A 50 µl reaction containing 11,000 cpm of a radiolabeled single-stranded DNA substrate and 10 µl of enzyme solution incubated for 4 hours at 37°C resulted in less than 5.0% release of TCA-soluble counts.

Double-Stranded Exonuclease Activity

A 50 µl reaction containing 5,000 cpm of a radiolabeled double-stranded DNA substrate and 10 µl of enzyme solution incubated for 4 hours at 37°C resulted in less than 0.5% release of TCA-soluble counts.

Endonuclease Activity

A 50 µl reaction containing 0.5 µg of pBR322 DNA and 10 µL of enzyme solution incubated for 4 hours at 37°C resulted in no visually discernible conversion to nicked circular DNA as determined by agarose gel electrophoresis.

E.coli 16S rDNA Contamination Test

Replicate 5 µl samples of enzyme solution were denatured and screened in a TaqMan qPCR assay for the presence of contaminating E.coli genomic DNA using oligonucleotide primers corresponding to the 16S rRNA locus. The acceptance criterion for the test is the threshold cycle count (Ct) produced by the average of 3 replicate no template control samples. Based on the correlation between the no template control Ct values, and standard curve data, the detection limit of this assay is <10 copies genome/sample.

QUALITY CONTROL ANALYSIS:

Unit Characterization Assay

Specific activity was measured using a 2-fold serial dilution method. Dilutions of enzyme were made in 1X reaction buffer. Reactions were incubated 10 minutes at 75°C, plunged on ice, and analyzed using the method of Sambrook and Russell.

SDS-Page (Physical Purity Assessment)

2.0 µl of concentrated enzyme solution was loaded on a denaturing 4-20% Tris-Glycine SDS-PAGE gel flanked by a broad-range MW marker and 2.0 µl of a 1:100 dilution of the sample. Following electrophoresis, the gel was stained and the samples compared to determine physical purity. The acceptance criteria for this test requires that the aggregate mass of contaminant bands in the concentrated sample do not exceed the mass of the protein of interest band in the dilute sample, confirming greater than 99% purity of the concentrated sample.

ORDERING INFORMATION

Product	Cat No.	Package
WizScript™ III Reverse Transcriptase (200U/µl)	W2303	10,000 U
	W2303-5	50,000 U

Technical Support



www.wizbiosolution.com
support@wizbiosolution.com
+82 70 7605 5066



Wizbiosolutions Inc.
A-802, Woolim Lions Valley 2, Sagimakgol-ro 45beon-gil 14,
Jungwon-gu, Seongnam-si 13209 Republic of Korea

FQP-806-02-114 (V.1.2)