

## DESCRIPTION

WizScript™ RT Master is a complete system for the efficient synthesis of first strand cDNA from RNA templates.

WizScript™ RT Master included MMLV RTase (RNase H-) which is an RNA-dependent DNA polymerase that is used in cDNA synthesis with long RNA templates. The lack of RNase H activity is important in this application in that RNase H activity will start to degrade template during long incubation times which are required for producing long cDNAs. RNase H minus RT enables preparation of long cDNAs and libraries containing a high percentage of full-length cDNA.

The kit is also supplied with both oligo dT and random primers. The oligo dT anneals selectively on the poly(A) tail of mRNA. Random primers do not require the presence of poly(A). Therefore, they can be used for transcription of the 5'-end regions of mRNA. Gene-specific primers may also be used with the kit. The first strand of cDNA can be directly used as a template in PCR.

## KIT CONTENTS

Contents	W2203
WizScript™ RT Master	1 ml
Oligo dT20 (50 pM)	100 µl
Random Hexamer (50 pM)	200 µl
RNase-Free Water	1 ml

- RT Master mix containing WizScript™ RTase, RNase inhibitor, stabilizer and reaction buffer containing optimized concentrations of MgCl<sub>2</sub>, dNTPs and DTT.

## APPLICATIONS

- First strand cDNA synthesis
- Full length cDNA library
- Real-time PCR
- PCR
- DNA labeling

## STORAGE CONDITIONS

Store all components at -20°C in a non-frost-free freezer.

## NOTE

Do not contaminate the WizScript™ RT Master with template RNA used in individual reactions.

## QUALITY CONTROL ANALYSIS:

In accordance with Wizbiosolutions Inc. ISO 13485-certified Quality Management System, each lot WizPure™ RT Master kit is tested against predetermined specifications to ensure consistent product quality.

## PROTOCOL

1. Prepare the following mixture in a microtube.

Component	Volume
RT Master mix	10 µl
Oligo dT (Random hexamer)	1 µl (2 µl)
Template RNA*	< 5 µg
RNase free water	up to 20 µl

- \* Notes : Recommended amounts of RNA template and primers for first-strand cDNA synthesis.

(1) RNAs : total RNA : 10 ng ~ 5 µg  
poly(A)+ RNA : 1 ng ~ 500 ng

(2) Primers : Oligo dT20 : 50 pM  
Random hexamer : 50 ~100 pM  
Gene-Specific Primer : 15 ~ 20 pM

2. Place tube in a thermal cycler programmed as follows:  
25°C / 10 min.  
42°C / 30 min.  
85°C / 5 min.  
4°C / Hold

3. Synthesized cDNA is immediately used as template for PCR or store at -20°C. The cDNA can be diluted with TE buffer and stored at -20°C.

## Notes

1. Isolation of poly(A)+ RNA from total RNA is not mandatory, however, doing so may improve the yield and purity of the final product.
2. RNA sample must be free of contaminating genomic DNA.
3. Unlike the oligo dT priming, which usually requires no optimization, the ratio of a random primer to RNA is critical in terms of the average length of cDNA synthesized in the reaction. Increasing the ratio of random primer/RNA will result in higher yield of shorter (~500 bp) cDNA, whereas decreasing this ratio will produce longer products.
4. The synthesized cDNA should be stored at -20°C.

## ORDERING INFORMATION

Product	Cat No.	Package
WizPure™ RT Master	W2203	100 rxn

## Technical Support



www.wizbiosolution.com  
support@wizbiosolution.com  
+82 70 7605 5066



**Wizbiosolutions Inc.**  
A-802, Woolim Lions Valley 2  
Sagimakgol-ro 45beon-gil 14  
Jungwon-gu, Seongnam-si  
13209 Republic of Korea