

## DESCRIPTION

WizPure™ HS-PCR 2X Master is ready-to-use Hot-start PCR pre-mixes are the innovation for convenience of your routine PCR.

The HS-PCR 2X Master contains an antibody-mediated hot-start Taq DNA Polymerase, MgCl<sub>2</sub>, dNTPs, enhancer and stabilizer.

The mixture is suitable for amplification of most of the DNA templates and highly processive 5'→3' DNA polymerase that lacks 3'→5' exonuclease activity and lacks a 3'→5' proofreading function. PCR reactions can be directly loaded onto an agarose gel without the additional need of loading buffer and dyes.

## KIT CONTENTS

| Contents                  | W1411 | W1411-5  |
|---------------------------|-------|----------|
| WizPure™ HS-PCR 2X Master | 1 ml  | 5 X 1 ml |

## APPLICATIONS

WizPure™ HS-PCR 2X Master is suitable and tested for amplification of genomic targets ranging from 100 bp to 4 kb and of episomal targets (lambda phage; plasmids) up to 10 kb under various reaction conditions.

- High through-put PCR
- Hot-start PCR
- Routine diagnostic PCR requiring high reproducibility
- DNA sequencing template preparation

## STORAGE CONDITIONS

Store all components at -20°C in a non-frost-free freezer.

## NOTE

Do not contaminate the WizPure™ HS-PCR 2X Master with primers and template DNA used in individual reactions. Thaw and mix all components thoroughly, spin down shortly and chill on ice.

## QUALITY CONTROL ANALYSIS:

In accordance with Wizbiosolutions Inc. ISO 13485-certified Quality Management System, each lot WizPure™ HS-PCR 2X Master kit is tested against predetermined specifications to ensure consistent product quality.

## PROTOCOL

This standard protocol applies to a single reaction where only template, primers, and water need to be added to the HS-PCR 2X Master mix. For multiple reactions, scale-up volume of reaction components proportionally. All reagents should be thawed on ice, gently mixed and briefly centrifuged before use.

1. Thaw reagents at room temperature. Mix thoroughly and then place on ice immediately after thawing.
2. Assemble reaction tubes on ice whenever possible to avoid premature, nonspecific polymerase activity.
3. The following table shows recommended component volumes:

| Component           | 20 µl reaction | Final Conc. |
|---------------------|----------------|-------------|
| HS-PCR 2X Master    | 10 µl          | 1X          |
| 10µM Forward Primer | 0.2 - 2.0 µl   | 0.1-1.0 µM  |
| 10µM Reverse Primer | 0.2 - 2.0 µl   | 0.1-1.0 µM  |
| Template DNA        | 1 - 5 µl       | < 250 ng    |
| Water, RNase-Free   | up to 20 µl    |             |

**NOTE:** In general, use greater than 0.5 µM primers for sensitivity and less than 0.5 µM for specificity.

**NOTE:** Recommended amount of template per PCR reaction:

- < 50 ng plasmid or
- < 500-1000ng genomic DNA or
- 2µl of a 100µl single plaque eluate or
- one single bacterial colony

4. Ensure reactions are mixed thoroughly by pipetting or gentle vortexing followed by a brief spin in a microcentrifuge.  
(Optional) Overlay reactions with one-half volume PCR-grade mineral oil when not using heated lid on thermal cycler.
5. Transfer tubes into a PCR instrument and run as following table.

| Step                 | Temp (°C) | Time         | Cycle   |
|----------------------|-----------|--------------|---------|
| Initial Denaturation | 95        | 5 min.       | 1       |
| Denature             | 95        | 10 - 60 sec. | 25 - 40 |
| Anneal               | 50 - 65   | 10 - 60 sec. |         |
| Extend               | 72        | 60 sec./kb   |         |
| Final Extension      | 72        | 5 min.       | 1       |

**NOTE:** Cycling conditions may need to be optimized, depending on different primer and template combinations. For example, raise the annealing temperature to prevent non-specific primer binding, increase extension time to generate longer PCR products.

6. After cycling, maintain the reactions at 4°C or store at -20°C until ready for analysis.

## ORDERING INFORMATION

| Product                   | Cat No. | Package |
|---------------------------|---------|---------|
| WizPure™ HS-PCR 2X Master | W1411   | 100 rxn |
|                           | W1411-5 | 500 rxn |

## Technical Support



[www.wizbiosolution.com](http://www.wizbiosolution.com)  
[support@wizbiosolution.com](mailto:support@wizbiosolution.com)  
 +82 70 7605 5066



**Wizbiosolutions Inc.**  
 A-802, Woolim Lions Valley 2  
 Sagimakgol-ro 45beon-gil 14  
 Jungwon-gu, Seongnam-si  
 13209 Republic of Korea